# CHROMABOND® HR-Xpert

the professional concept of innovative SPE phases

CHROMABOND® HR. Xperl

available
in
ana 85 mm

polymer-based RP-phases and mixed-mode ion exchange phases

**MACHEREY-NAGEL** 

www.mn-net.com





### CHROMABOND® HR-Xpert

... the innovative concept of five polymer-based RP- and mixed-mode ion exchange phases for SPE:

> CHROMABOND® HR-X

> CHROMABOND® HR-XC

> CHROMABOND® HR-XA

> CHROMABOND® HR-XCW

> CHROMABOND® HR-XAW


hydrophobic PS/DVB copolymer strong mixed-mode cation exchanger strong mixed-mode anion exchanger weak mixed-mode cation exchanger weak mixed-mode anion exchanger

#### ... these innovative SPE phases offer:

state-of-the-art spherical polymer

- $\Rightarrow$  different particle sizes (45  $\mu$ m and 85  $\mu$ m) suitable depending on the sample volume and matrix
- \Rightarrow broad spectrum of applications with special suitability for enrichment of pharmaceuticals from biological matrices
- ideal flow properties due to low content of particulate matter
- optimized pore structure and high specific surface
  - high loadability and outstanding elution properties
  - ⇒ low solvent consumption
  - rapid, economical analyses
- high-purity adsorber material
  - allows highest reproducibility with extremely low blind values
  - > reliable analyses at ultra trace level
  - no method adaptation for new batches necessary



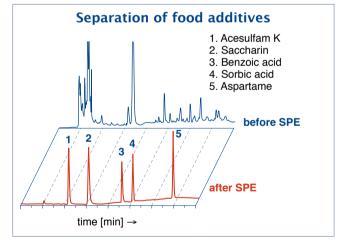

#### In modern analyses SPE fulfills 3 important tasks:

- => enrichment of analytes up to factor 10.000 - increase of chromatographic sensibility/improved limits of detection
- > removal of interfering compounds protection of subsequent analyses like HPLC, GC, TLC, MS, UV or IR spectroscopy, ...
- changing of an analyte's matrix matrix which is simpler or more suitable for subsequent analyses


# ... the HR-X*pert* concept guarantees:

- RP- and mixed-mode SPE-phases with distinct ion exchange and reversed-phase properties your benefit ⇒ excellent enrichment of neutral, acidic and basic compounds
- omodern, spherical support-polymer with optimized pore structure and high surface your benefit ⇒ good reproducibility, reliable and cost-efficient analysis
- possibility for more aggressive washing procedures for matrix removal your benefit ⇒ cleaner samples and protection of your HPLC and GC-machines
- quantification of analytes also from heavily contaminated samples your benefit ⇒ lower limits of detection also for critical matrices

CHROMABOND® HR-Xpert is the perfect combination for all tasks in sample preparation




### **Chemical structures**

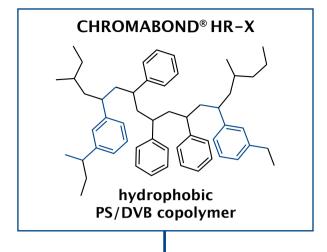

#### CHROMABOND® HR-XCW


weak acidic cation exchanger

#### CHROMABOND® HR-XA

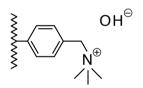


strong basic anion exchanger

#### CHROMABOND® HR-XAW

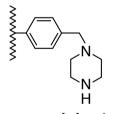


weak basic anion exchanger

#### CHROMABOND® HR-XC

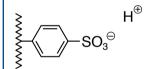

strong acidic cation exchanger

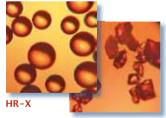

conventional PS/DVB copolymer

HR-X: spherical base material for efficient enrichment and ideal flow behavior

### Similary phases:

**CHROMABOND® HR-X:** Oasis® HLB, Strata™ X, Nexus, ENVI-Chrom P

CHROMABOND® HR-XC: Oasis® MCX, Strata™ X-C, StyreScreen® DBX, HyperSep™ Retain™-CX
CHROMABOND® HR-XA: Oasis® MAX, Strata™ X-A, HyperSep™ Retain™-AX, StyreScreen® QAX

CHROMABOND® HR-XCW: Oasis® WCX, Strata™-X-CW
CHROMABOND® HR-XAW: Oasis® WAX, Strata™-X-AW
(Trademarks of Waters, Phenomenex, Varian, Thermo Fisher Scientific Inc. and UCT)


www.mn-net.com \_\_\_\_\_\_ 3 \_\_\_\_

MN




# 2 particle sizes - 1 goal ... with HR-X pert to the ideal application

for different applicative requirements the particle sizes complement each other perfectly

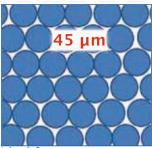

#### ideal for:

- smaller sample volumes
- smaller amount of sorbent
- ⊕ lower elution volumes

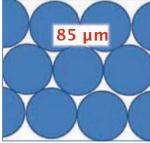

- voluminous or viscous samples/ strong matrix load
- operation without vacuum (e.g. for volatile analytes)
- higher amount of sorbent without pressure rise

#### Comparison of 45 μm and 85 μm

Larger number of particles per volume at 45 µm by about half radius

ightfold particle number at same amount of sorbent

Specific surface is equal for both particle sizes


🕏 freely accessible external surface is at 45 μm particles much larger


More compact sorbent package: Enhanced interaction of the analytes with the sorbent

improved extraction results

#### Ideal elution characteristics

Method: 1 mL column with 30 mg CHROMABOND® HR-X, 1 mL standard solution (1 mg/mL hexobarbital); drying; elution in portions of 100 µL with methanol, according to MN Appl. No. 305490

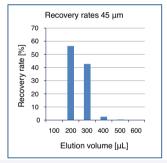

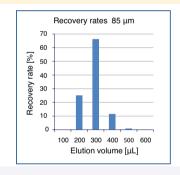

at 45 μm · faster elution · lower elution volumes needed

### Breakthrough behavior in enrichment

Method: 1 mL column with 15 mg CHROMABOND® HR-X; apply portions of 1 mL standard solution (250 µg/mL hexobarbital in water); collect the eluate, according to MN Appl. No. 305480

#### **45 μm** (red):


Analyte is completely retained up to the fraction 10.

#### 85 µm (blue):

Small amounts even break through at fraction 4.

### Comparison of 45 μm and 85 μm

45 µm are smaller particles and provide better enrichment of sorbent and a superior breakthrough behavior.

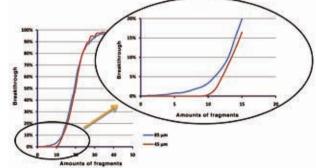


O When using larger amounts of sorbent this effect is less distinct. Also using 85 μm particles sufficient contact of analyte and sorbent is given. So, 45 µm particles are ideal for small sample and elution volumes, while for large amounts of sample or sorbent 85 µm particles provide advantages due to with better flow properties.



# 3 paths - 1 goal ... cleaner samples by HR-Xpert:

Depending on the character of the analytes HR-Xpert offers the suitable adsorbent and optimal method for sample preparation, cleaning and concentration.

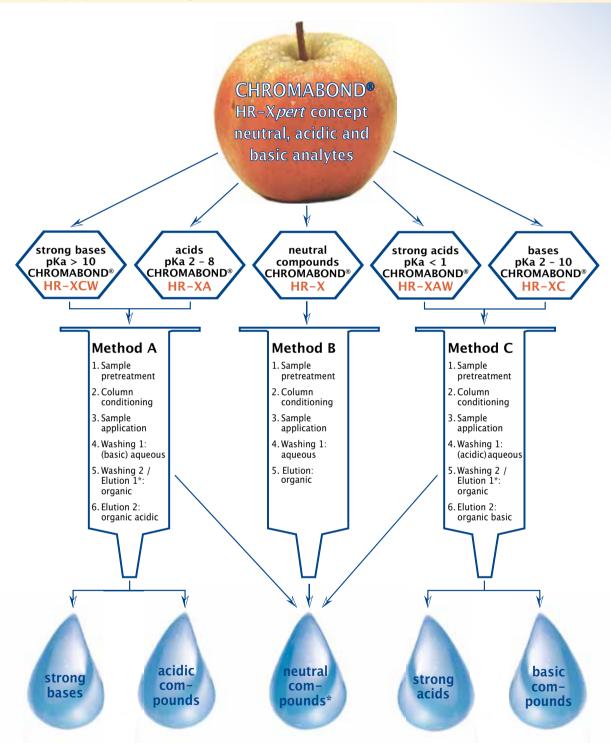

#### \* under organic washing and elution conditions the following compounds will be also eluted:

HR-X: polar compounds as organic acids and bases

HR-XC/HR- XCW: acidic compounds/impurities HR-XA/HR- XAW: basic compounds/impurities

www.mn-net.com

(MN)



# HR-X

# spherical, hydrophobic polystyrene-divinylbenzene adsorbent resin

- hydrophobic polystyrene-divinylbenzene copolymer pH stability 1-14
- high-purity material with highest reproducibility and lowest blank values due to a novel manufacturing process
- spherical particles, size 45 μm and 85 μm (standard);
   pore size 55-60 Å, very high surface 1000 m²/g
   capacity 390 mg/g (caffeine in water)
- excellent recovery rates especially for the enrichment of pharmaceuticals / active ingredients due to the spherical structure of the particles, very homogeneous surface, and optimized pore structure
- recommended application:
  pharmaceuticals / active inquired.

pharmaceuticals / active ingredients from tablets, creams and water / waste water


drugs and pharmaceuticals from urine, blood, serum and plasma trace analysis of pesticides, herbicides, phenols, PAHs and PCBs from water

# Standard protocol for CHROMABOND® HR-X

MN Appl. No. 304310

Column type: CHROMABOND® HR-X/3 mL/200 mg, REF 730931

Sample pretreatment: individual sample preparation with reference to analytes and matrix

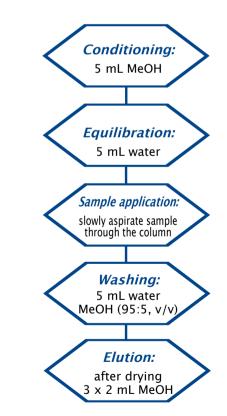

Further analysis: if necessary, evaporate and redissolve in a suitable solvent; HPLC or GC

These conditions are a starting point for SPE method development. Further optimisation may be required to improve results.

#### Sulfonamides from serum

MN Appl. No. 304220

Column type: CHROMABOND® HR-X/3 mL/200 mg

REF 730931

Sample: 2 µg/ml each in serum

Column conditioning: 5 mL MeOH, 5 mL dest. water

Sample application: 1 mL spiked serum

Column washing: 5 mL water - MeOH (95:5, v/v)

Elution: after drying 3 x 2 mL MeOH

#### Recovery rate [%]:

Compound	HR-X	Oasis® HLB	Strata™ X
Sulfanilamide	66	62	63
Sulfadiazine	107	101	108
Sulfamerazine	114	111	111
Sulfadimidine	94	86	89
Succinylsulfathiazole	70	43	48

#### Further analysis:

HPLC see MN Appl. No. 117880

Column: 125/4 mm

NUCLEODUR® C<sub>18</sub> Gravity, 5 µm

Eluent: MeOH - 0.1 % TFA (20:80, v/v)


Flowrate: 1.0 mL/min

Temperature: 22 °C

Detection: UV, 230 nm

#### Peaks:

- 1. Sulfanilamide
- 2. Sulfadiazine
- 3. Sulfamerazine
- 4. Sulfadimidine
- 5. Succinylsulfathiazole

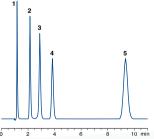




MN



### **Highest reproducibility**

- √ within each batch
- √ from batch to batch

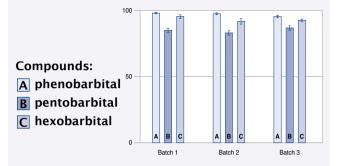

#### **Barbiturates from serum**

MN Appl. No. 304290

Column type: CHROMABOND® HR-X/3 mL/200 mg REF 730931

Sample: 100 ng/ml each in serum

Column conditioning: 5 mL MeOH, 5 mL dist. water

Sample application: 1 mL spiked Serum

Column washing: 5 mL water

Elution: after drying 3 x 2 mL MeOH

Further analysis: HPLC on NUCLEODUR® 100-5 C, ec,

see MN Appl. No.117820

#### Pesticides from water

MN Appl. No. 304250 and 304260 GIU GmbH, Teningen

Column type: CHROMABOND® HR-X/3 mL/200 mg

REF 730931

Sample: 500 ng/L each in water; pH value adjusted Column conditioning: 10 mL MeOH, 10 mL water Sample application: slowly aspirate 1 L sample

through the column Drying: stream of nitrogen

Elution: 5 mL MeOH - THF (1:1, v/v)

#### Recovery rate [%]:

Compound	pH 7	Compound	pH 2
Desisopropylatrazine	90	Metamitrone	86
2,4-Dichlorbenzamide	95	Quinmerac	90
Desethylatrazine	89	Chloridazone	93
Hexazinone	95	Picloram	83
Bromacil	103	Metribuzine	84
Simazine	91	Cyanazine	83
Desethylterbutylazine	89	Metabenzthiazuron	94
Atrazine	88	Chlortoluron	91
Metalaxyl	97	Isoproturon	89
Metazachlor	93	Diuron	91
Propazine	88	Dimethenamid-P	89
Terbuthylazine	86	Linuron	94
Metolachlor	97	Epoxiconazole	86
		Penconazole	90
		Alachlor	93
		Propiconazole-1	89
Further analysis:		Flufenacet	91

HPLC e.g. on NUCLEODUR® 100-3 C. ec see MN Appl. No. 118610 or 120485

# **Drugs from water**

MN Appl. No. 304240

Column type: CHROMABOND® HR-X/3 mL/200 mg

REF 730931

Sample: each 1 µg/L in water

Column conditioning: 5 mL MeOH, 5 mL water Sample application: slowly aspirate 500 mL water

(pH 3) through the column Column washing: 5 mL water

*Elution:* after drying 3 x 2 mL acetonitrile

#### Pacovary rata [%]

necovery rate [/0].		
Compound	HR-X	Strata™ X
Ketoprofen	98	92
Ibuprofen	91	93
Pentobarbital	99	95
Meclofenamic acid	92	93
Protriptyline	63	45
Nortriptyline	53	39

Further analysis: HPLC see MN Appl. No. 121690

Column: 125/4 mm

NUCLEODUR® C<sub>18</sub> Gravity, 5 µm

Eluent: A) 20 mmol KH<sub>2</sub>PO<sub>4</sub>, pH 7

B) acetonitrile

Gradient:  $20 \% B \rightarrow 50 \% B in 25 min$ 

Flow rate: 0.5 mL/min

Temperature: 22 °C


Detection: UV, 215 nm

#### Peaks:

1. Ketoprofen 2. Ibuprofen

3. Pentobarbital

6. Nortriptyline

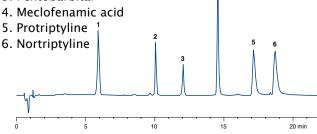

www.mn-net.com



# HR-XC

- strong acidic benzenesulphonic acid cation exchanger exchange capacity 1.0 meg/g, pKa <1
- base material polystyrene-divinylbenzene copolymer pH stability 1-14
- high purity material, highest reproducibility and lowest blank values due to an optimized production process
- spherical particles, size 45 μm and 85 μm (standard); pore size 65-75 Å, very large specific surface 800 m<sup>2</sup>/g; pore volume 1.4 cm<sup>3</sup>/g, RP capacity 300 mg/g (caffeine in water)
- outstanding recovery rates especially for the enrichment of basic analytes

#### strong cation exchanger

recommended application: basic active ingredients from heavily matrix-contaminated samples like e.g. urine, plasma, fungicides from food melamine from milk products basic analytes like e.g. amines

#### Standard protocol for CHROMABOND® HR-XC

MN Appl. No. 304790

Column type: CHROMABOND® HR-XC/3 mL, 200 mg, REF 730952

Sample pretreatment: individual sample preparation with reference to analytes and matrix

> Conditionina: 5 mL MeOH **Equilibration:**

5 mL water

Sample application: slowly aspirate sample through the column

Washing1: 2 mL 0.1 mol/L HCl in water

Washing 2:/Elution1: 2 mL MeOH (elution of neutral and acidic compounds)

Elution2: after drying 5 mL MeOH/5 % NH, (elution of basic compounds)

Further analysis: if necessary, evaporate and redissolve in a suitable solvent; HPLC or GC

These conditions are a starting point for SPE method development. Further optimisation may be required to improve results.

### Fractionation of acidic, neutral and basic analytes from serum

bases with pKa 2 - 10

MN Appl. No. 304780

Column type: CHROMABOND® HR-XC/3 mL, 200 mg, REF 730952

Sample: 1 mL spiked matrix, acidified with 200 µL 2 % H<sub>3</sub>PO<sub>4</sub>

Column conditioning: 5 mL MeOH, then 5 mL water


Sample application: 1.2 mL spiked serum

Washing: 2 mL 0.1 N HCl in water

Elution: a) 2.5 mL MeOH

Fraction A: neutral and acidic analytes

b) 5 mL MeOH/10 % NH, Fraction B: basic analytes

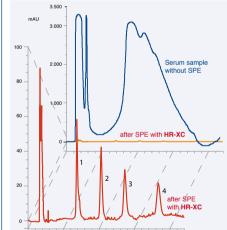

#### Peaks:

- 1. Doxepin
- 2. Imipramine
- 3. Amitriptyline
- 4. Trimipramine

#### Recovery rates [%]:

Fraction A: acidic Fraction B: basic analytes and neutral analytes

Co	mpound	HR-XC	Compound	HR-XC	Oasis® MCX	Strata™ X-C
Sup	orofen	108	Doxepin	101	68	82
Naj	oroxen	85	Imipramine	95	71	85
Tol	metin	73	Amitriptyline	94	72	78
Phe	nobarbital	108	Trimipramine	92	70	81
Ind	omethacin	33				_
He	xobarbital	80				

Further analysis: HPLC on NUCLEODUR® C18 e.g. C8 Gravity; see MN Appl. No. 122230 and 118520

(MN



# HR-XA

- strong basic quaternary ammonium anion exchanger
- exchange capacity 0.25 meq/g, pKa ~ 18

  base material polystyrene-divinylbenzene copolymer
  pH stability 1 14
- high purity material with highest reproducibility and lowest blank values due to an optimized production process
- spherical particles, size 45 μm and 85 μm (standard);
   pore size 55 65 Å, very large specific surface 850 m²/g; pore volume 1.4 cm³/g, RP capacity 350 mg/g (caffeine in water)
- outstanding recovery rates especially for the enrichment of acidic analytes

#### strong anion exchanger

recommended application:
acidic active ingredients from
heavily matrix-contaminated
samples like e. g. urine, plasma,
serum
phenolic acids
acidic herbicides
weak/medium-strength acids with
pKa 2 - 8



# Acidic Pharmaceuticals from serum

MN Appl. No. 305000

Column type: CHROMABOND® HR-XA/3 mL,

200 mg, REF 730951

Sample: 5 mL spiked serum (1 µg/ml each), adjusted

basic with 1 N NaOH

Column conditioning: 5 mL MeOH

Equilibration: 5 mL water

Sample application: apply sample

Washing: a) 2.5 mL water

b) 2.5 mL 0,1 N NaOH

c) 5 mL MeOH

Elution: after drying, 5 mL MeOH /1 % formic acid

Further analysis: HPLC on

NUCLEODUR® C18 Gravity; see MN Appl.-No. 122840

Recovery rate [%]:

HR-XA	Oasis® MAX
90	85
104	123
98	69
106	98
88	58
69	89
95	94
92	93
	90 104 98 106 88 69

# Further applications on CHROMABOND® HR-XA

- Appl. No. 304980 Acidic Herbicides
- Appl. No. 305020 Fractionation of Acidic and Basic Analytes from serum
- Appl. No. 304990 Phenolic Acids


www.mn-net.com/apps

# Standard protocol for CHROMABOND® HR-XA

MN Appl. No. 304970

Column type: CHROMABOND® HR-XA/3 mL, 200 mg, REF 730951

Sample pretreatment: individual sample preparation with reference to analytes and matrix

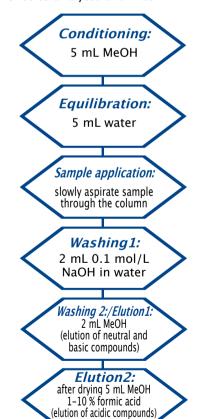

Further analysis: if necessary, evaporate and redissolve in a suitable solvent: HPLC or GC

These conditions are a starting point for SPE method development. Further optimisation may be required to improve results.

-(MN



#### HR-XCW

- weak acidic carboxylic acid cation exchanger exchange capacity >0.7 meg/g, pKa ~ 5
- base material spherical PS/DVB copolymer pH stability 1-14
- high purity material, highest reproducibility and lowest blank values due to an optimized production process
- spherical particles, size 45 μm and 85 μm (standard); pore size 50-60 Å very large specific surface 850 m²/g; pore volume 1.2-1.4 cm³/g RP capacity 350 mg/g (caffeine in water)
- outstanding recovery rates especially for enrichment of strongly basic analytes

#### weak cation exchanger

recommended application:

basic compounds like quaternary amines active ingredients from heavily matrix-contaminated samples like e.g. urine, plasma, serum strong bases with pKa > 10

# Standard protocol for CHROMABOND® HR-XCW

MN Appl. No. 305300

Column type: CHROMABOND® HR-XCW/3 mL, 200 mg, REF 730739

Sample pretreatment: individual sample preparation with reference to analytes and matrix

#### Conditioning: 5 mL MeOH

#### **Equilibration:**

5 mL water

#### Sample application:

slowly aspirate sample through the column

#### Washing1:

2 mL acidified water (1-5 % formic acid)

### Washing 2:/Elution1:

2 mL MeOH (elution of neutral and acidic compounds)

#### Elution2:

after drying 2 x 2 mL MeOH / 1-5 % formic acid (elution of strongly basic compounds)

Alternatively elution 2 can be achieved by basic methanol ( $\mathrm{NH_3}$ ), e.g. for primary to tertiary amines. Deprotonation of the analytes interrupts the interaction with the cation exchange resin.

Further analysis: if necessary, evaporate and redissolve in a suitable solvent; HPLC or GC

These conditions are a starting point for SPE method development. Further optimisation may be required to improve results.

### **Tricyclic Antidepressants**

MN Appl. No. 305340

Column type: CHROMABOND® HR-XCW/3 mL, 60 mg REF 730735

Sample: 250  $\mu$ L spiked Serum, diluted with 1 mL 10 %

formic acid in water

Conditioning: 3 mL MeOH Equilibration: 3 mL water

Sample application: slowly aspirate sample through

the column

Washing: 1 mL 5 % formic acid in water, then

1 mL MeOH

*Elution:* after drying by vaccum (15 min)

3 mL 5 % formic acid in MeOH

Further analysis:

evaporate and redissolve in a suitable solvent for HPLC on NUCLEODUR® C8; Gravity see

MN Appl. No. 118520

Recovery rate [%]:

#### Compound HR-XCW HR-XC\* PCA\*\* Oasis® WCX Doxepine 79 5 11 41 79 9 67 **Imipramine** 20 Amitriptyline 91 9 14 46 **Trimipramine** 98 7 27

\* HR-XC: Basic analytes can not be eluted with slightly acidic organic conditions from the strong cation exchanger CHROMABOND® HR-XC, because the eluting power is not sufficient to dissociate the interaction with the ion exchanger. However, with the usage of basic methanol a complete elution can be achieved (please see also MN Appl. No. 304780).

\*\* PCA: Due to the missing RP interactions of silica based weak cation exchanger, CHROMABOND® PCA gives only a small enrichment elution of the analytes.





#### HR-XAW

- weak basic secondary and tertiary ammonium anion exchanger exchange capacity > 0.5 meg/g, pKa ~ 6
- base material spherical PS/DVB copolymer pH stability 1-14
- high purity material with highest reproducibility and lowest blank values due to an optimized production process
- spherical particles, size 45 µm and 85 µm (standard); pore size 55-65 Å very large specific surface 850 m<sup>2</sup>/g; pore volume 1.2-1.4 cm<sup>3</sup>/g RP capacity 350 mg/g (caffeine in water)
- outstanding recovery rates especially for enrichment of acidic analytes

#### weak anion exchanger

recommended application: perfluorinated surfactants acidic compounds like sulfonates active ingredients from heavily matrix-contaminated samples like e.g. urine, plasma, serum strong acids with pKa < 1



Impregnated with fluorosurfactents?

# Surfactants from water MN Appl. No. 305140

Column type: CHROMABOND® HR-XAW/3 mL, 60 mg, REF 730747

Sample: 500 mL water, spiked with 1 mL standard solution (20 µg/L of each compound)

Conditioning: 2 mL MeOH + 5 % ammonia, then

2 mL methanol Equilibration: 2 mL water

Sample application: slowly aspirate sample through

the column

Washing: 1. 2 mL water

2. 2 mL acetone - acetonitrile formic acid (50:50:1, v/v/v),

3. 2 mL MeOH

Elution: after drying by vaccum 2 mL methanol + 5 % ammonia

#### Futher analysis:

evaporate to dryness in a stream of nitrogen under slight heating, and redissolve in a suitable solvent for HPLC on NUCLEODUR® Sphinx RP; see MN Appl. No. 123340

#### Recovery rate [%]:

recovery rate [/o].		
Compound	HR-XAW	Oasis® WAX
Perfluoropropionic acid (PFPrA)	103	76
Perfluoropentanoid acid (PFPeA)	94	81
Perfluorohexanoic acid (PFHxA)	94	91
Perfluoroctanoic acid (PFOA)	95	95
Perfluoroctane sulfonate K salt (PFOS)	81	80
Perfluorododecanoic acid (PFDoDA)	82	34

Application in accordance with **DIN 38407-42.** 

#### Standard protocol for CHROMABOND® HR-XAW

MN Appl. No. 305200

Column type: CHROMABOND® HR-XAW/3 mL, 200 mg, REF 730748

Sample pretreatment: individual sample preparation with reference to analytes and matrix



Equilibration: 5 mL water

Sample application:

slowly aspirate sample through the column

### Washing 1:

25 mM ammonium acetate in water

Washing 2:/Elution1: 2 mL MeOH (elution of neutral and basic compounds)

Elution2: after drying 2 x 2 mL MeOH 1-5 % NH<sub>3</sub> (elution of strongly

acidic compounds)

Alternatively elution 2 can be achieved by acidified methanol (formic acid). Protonation of the analytes interrupts the interaction with the anion exchange resin.

Further analysis: if necessary, evaporate and redissolve in a suitable solvent; HPLC or GC

These conditions are a starting point for SPE method development. Further optimisation may be required to improve results.

www.mn-net.com 11-

**《MN**)



#### Alkaloids from serum

MN Appl. No. 304300

Column type: CHROMABOND® HR-X/3 mL/200 mg

REF 730931

Sample pretreatment: Serum samples are spiked with alkaloids (each 2 µg/mL).

Conditioning: 5 mL MeOH

Equilibriation: 5 mL water (Do not let run the column

dry!)

Sample application: 1 mL spiked serum sample is

passed through the column by vacuum

Washing: 5 mL water Drying: with nitrogen or air

Elution: after drying 3 x 2 mL acetone

Further analysis: HPLC with NUCLEODUR  $^{\rm @}$  C  $_{\rm 18}$  Gravity, according to MN Appl. No. 117950

**Recovery rates [%]:** 

Compound	HR-X
Atropine	99
Papvarine	97
Noscapine	95
Strychnine	94
Quinine	60

#### Fungicides in fruit juice

MN Appl. No. 304740

Column type: CHROMABOND® HR-XC/3 mL/200 mg

REF 730952

Sample pretreatment: 5 mL spiked apple juice

 $(c=1 \mu g/mL) + 0.5 mL 0.1 N NaOH$ 

Conditioning: 5 mL MeOH

Equilibriation: 5 mL 2 % NH<sub>3</sub> (Do not let run the

column dry!)

Sample application: prepared sample is passed

through the column by vacuum.

Washing:

a) 2 mL 2 % NH3

b) 2 mL 30 % MeOH/5 % NH<sub>3</sub>

c) 2 mL 0.1 N HCl

d) 2 mL MeOH

Drying: with nitrogen or air Elution: 4 mL MeOH/5 % NH,

Evaporation to dryness and reconstitution with 1 mL

of mobile phase from subsequent HPLC

Further analysis: HPLC with NUCLEODUR® C1. Gravity,

according to MN Appl. No. 122200

**Recovery rates [%]:** 

Compound	HR-XC
Carbendazime	89
Thiabendazole	92
<u> </u>	

# **Extraction of melamine from**

MN Appl. No. 304920

Column type: CHROMABOND® HR-XC/3 mL/200 mg

REF 730952

Step 1: Add 1 mL 1N HCl to 10 mL spiked milk  $(1 \mu g/mL \text{ or } 0.1 \mu g/mL)$ ; then add 10 mL methylene chloride.

Step 2: vortex to mix, centrifuge for 15 min at 4500 rpm, remove the aqueous layer add 5 mL 0.1 N HCl to the organic layer

repeat twice step 2

At least combine all aqueous phases for SPE.

Conditioning: 5 mL MeOH

Equilibriation: 5 mL water (Do not let run the column

dry!)

Sample application: The prepared sample is passed through the column by gravity or low vacuum.

Washing: 5 mL 0,1 N HCl + 5 mL MeOH

Drying: 10 min vacuum

Elution: 2 x 2.5 mL MeOH/5 % NH<sub>3</sub>

Evaporation with nitrogen stream and reconstitution with a suitable solvent for the subsequent analysis Further analysis: HPLC with NUCLEODUR® 100-5 C. ec, according to MN Appl. No. 122670

Recovery rates [%]:

Compound	HR-XC
1 µg melamine in 1 mL milk	99
0.1 µg melamin in 1 mL milk	88

#### **Extraction of** campher sulfonic acid from serum

MN Appl. No. 305320

Column type: CHROMABOND® HR-XAW/3 mL/60 mg REF 730747

Sample pretreatment: Stock solution: Camphor sulfonic acid in water (c=0.5 mg/mL)

Standard solution: 5 µL stock solution is diluted to  $10 \text{ mL} (c=0.25 \mu g/mL)$ 

Sample solution: 5 mL serum is spiked with 5 µL stock solution and diluted with 4 % H<sub>3</sub>PO<sub>4</sub> to 10 mL

Conditioning: 2 mL MeOH

Equilibriation: 2 mL water (Do not let run the column

dry!)

Sample application: 2 mL sample solution is passed through the column by vacuum.

Washing: 2 mL ormic acid in water, then 2 mL MeOH

**Drying:** with vacuum

Elution: 2 mL MeOH + 5 % ammonia

Evaporation to dryness (mild nitrogen stream, slightly warming), reconstitution in suitable eluent

Further analysis: HPLC-MS mit NUCLEODUR® C<sub>18</sub>

Gravity, according to MN Appl. No. 123520 Recovery rate: 98.9 %



#### Allantoin in cosmetic products

MN Appl. No. 305440

Column type: CHROMABOND® HR-XA/3 mL/60 mg

REF 730950

Sample pretreatment: mix 1 q of cosmetic product with 100 mL of water and adjust pH to 10 - 11 with conc. ammonia solution.

Conditioning: 1 mL MeOH

Equilibriation: 1 mL ammonium hydroxide solution

 $(w(NH_3) = 5 \%)$ 

Sample application: slowly force or aspirate 1 mL

sample through the column.

Washing: 1 mL ammonium hydroxide solution  $(w(NH_3) = 5 \%$ , then 1 mL MeOH (do not dry the col-


umn completely!)

Elution: 2 x 600 µL HCI (w(HCI) = 0.6 %)

Fill up to 10 mL with acetonitrile/30 mM ammonium chloride, pH 3.0 (80:8, v/v) for HPLC analysis.

Further analysis: HPLC with NUCLEODUR® HILIC,

according to MN Appl. No. 305446

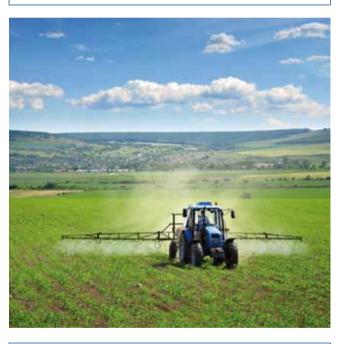

#### **Desalting and selective** Elution of peptides and proteins

MN Appl. No. 305290

Column type: CHROMABOND® HR-X/1 mL/30 mg REF 730934

Conditioning: 1 mL acetonitrile

Equilibriation: 1 mL de-ionized water (0.1 % TFA)

Sample application: Slowly force or aspirate the sample

through the column.

Washing: 2 x 1 mL de-ionized water (0.1 % TFA)

Elution: 250 µL of

40:60 ACN/water (acidic peptides) 70:30 ACN/water (neutral peptides)

70:30 ACN/1.0 % formic acid (basic peptides)

#### Acidic herbicides

MN Appl. No. 304980

Column type: CHROMABOND® HR-XA/3 mL/200 mg

REF 730951

Sample pretreatment: 1 µg/mL sample adjusted on

basic pH with 1 N NaOH Conditioning: 5 mL MeOH

Equilibriation: 5 mL water (Do not let run the column

dry!)

Sample application: prepared sample is passed

through the column by vacuum.

a) 2.5 mL 50 mM NaOAc

b) 2.5 mL MeOH

Drying: with nitrogen or air

Elution: 5 mL MeOH/10 % formic acid

Evaporation to dryness and reconstitution with 1 mL

of mobile phase from subsequent HPLC

Further analysis: HPLC with NUCLEODUR®  $\rm C_{18}$  Gravity, according to MN Appl. No. 122820

**Recovery rates [%]:** 

Compound	HR-XA	Oasis® WAX
Dicamba	87	78
Bentazon	79	72
2,4-D	82	80
MCPA	76	72

#### Extraction of paraquat and diquat from water

MN Appl. No. 305370

Column type: CHROMABOND® HR-XCW/3 mL/60 mg

REF 730735

Conditioning: 1 mL MeOH

Equilibriation: 1 mL water (Do not let run the column

dry!)

Sample application: 1 mL spiked water (0.5 µg/mL)

Washing: 1 mL water Drying: with nitrogen or air

Elution: 1 mL acetonitrile/water + 2 % formic acid

(1:1, v/v)

Evaporation and reconstitution is suitable solvent Further analysis: HPLC with NUCLEODUR® HILIC,

according to MN Appl. No. 123060

Recovery rates [%]:

Compound	HR-XA
Paraquat	103
Diquat	113

www.mn-net.com



# CHROMABOND® HR-X

### hydrophobic RP-phase

					,		P	
	Volume		Ads	orbent weig	ht			Pack of
	CHROMABOND® HR-X polypropylene columns (85 μm)							
		30 mg	60 mg	100 mg	200 mg	500 mg	1 g	
	1 mL 3 mL 6 mL 15 mL	730934	730936	730935	730931 730938	730937 730939 730940	730941	30 30 30 20
	<b>CHROM</b>	ABOND® HR-X	oolypropylene co	lumns (85	μm) · BIGpa	cks		
	3 mL 6 mL				730931.250 730938.250	730939.250	)	250 250
	<b>CHROM</b>	ABOND® HR-X I	oolypropylene co	lumns (45	μm)· NEW			
	1 mL 3 mL	730934P45	7 730936P45	'30935P45	730931P45			30 30
	CHROM	1ABOND® LV-HR	-X (85 μm)					
		30 mg	60 mg		200 mg			
	15 mL	732130	732131		732132			30
	<b>CHROM</b>	MABOND® MULTI	96 HR-X					
		96 x 10 mg (45 μm)	96 x 25 mg (45 µm)		50 mg μm)	96 x 10 (85 μ		
		738530.010M	738530.025M	73853	0.050M	738530.	100M	1
	<b>CHROM</b>	MABOND® HR-X a	adsorbent (85 µn	1)				
Charles and the control of the contr						7306	63	20 g

# CHROMABOND® HR-XC

## strong cation exchanger

	Volume			Adsorben	t weight			Pack of
	CHROMABOND® HR-XC polypropylene columns (85 µm)							
		30 mg	60 mg	100 mg	150 mg	200 mg	500 mg	
	1 mL	730969		730049				30
	3 mL		730956			730952	730953	30
	6 mL				730957		730955	30
CHROMABOND® HR-XC polypropylene columns (45 μm) · NEW								
	1 mL	730969P45		730049P45	-			30
	3 mL		730956P45			730952P45		30
	CHROMAFIX® HR-XC cartridges (85 µm)							
		Size	S		M		L	
	adsorb	ent weight Ø	155 mg		240 mg		500 mg	
			731755		731756		731757	50
70 PM	CHRON	MABOND® HR	-XC adsorbe	ent (85 µm)				
9343333333355 944333333335 9443333333335 9443333333333						730	0664	100 g

# CHROMABOND® HR-XA

# strong anion exchanger

								_
Volume			Adsorbent weight					Pack of
	CHROMABOND® HR-XA polypropylene columns (85 μm)							
		30 mg	60 mg	100 mg	150 mg	200 mg	500 mg	
	1 mL	730968		730727				30
	3 mL		730950			730951	730954	30
	6 mL				730958		730966	30
	CHROMABOND® HR-XA polypropylene columns (45 μm) · NEW							
	1 mL	730968P45		730727P45				30
	3 mL		730950P45			730951P45		30
	CHROMAFIX® HR-XA cartridges (85 μm)							
		Size	S		M		L	
W	adsorb	ent weight Ø	155 mg		240 mg		500 mg	
			731768		731769		731770	50
-088883bb	CHRON	MABOND® HR	-XA adsorbe	nt (85 µm)				
633333333333355						7306	571	100 g



# CHROMABOND® HR-XCW

# weak cation exchanger

Volume			Adsorbent weight					Pack of
	CHROMABOND® HR-XCW polypropylene columns (85 μm)							
		30 mg	60 mg	100 mg	150 mg	200 mg	500 mg	
	1 mL	730731		730733				30
T	3 mL		730735			730739	730741	30
	6 mL				730737		730743	30
CHROMABOND® HR-XCW polypropylene columns (45 µm) · N								
	1 mL	730731P45		730733P45				30
	3 mL		730735P45			730739P45		30
	CHROMAFIX® HR-XCW cartridges (85 μm)							
	-	Size nt weight Ø	<b>S</b> 155 mg		<b>M</b> 240 mg		<b>L</b> 500 mg	
			731774		731775		731776	50
	CHROMA	ABOND® HR-	XCW adsor	bent (85 µm)	)			
							730674	100 g

# CHROMABOND® HR-XAW

# weak anion exchanger

Volume			Adsorbent weight					Pack of
	CHROMABOND® HR-XAW polypropylene columns (85 μm)							
		30 mg	60 mg	100 mg	150 mg	200 mg	500 mg	
	1 mL	730728		730729				30
T	3 mL		730747			730748	730744	30
	6 mL				730749		730745	30
	<b>CHROM</b>	ABOND® HR-	-XAW polyp	ropylene col	umns (45	μm) · NEW		
	1 mL	730728P45		730729P45				30
	3 mL		730747P45			730748P45		30
	CHROMAFIX® HR-XAW cartridges (85 μm)							
	9	Size	S		M		L	
	adsorbe	nt weight $arnothing$	155 mg		240 mg		500 mg	
			731771		731772		731773	50
400000000000000000000000000000000000000	CHROMA	ABOND® HR-	-XAW adsor	rbent (85 µm	)			
(1) (1) (1) (1) (1) (1) (1) (1) (1) (1)							730673	100 g

# CHROMABOND® HR-Xpert

# development kits

			Pack of
CHROMABOND® polypropylene columns			
10 x HR-X and	Kit I 3mL/60mg	Kit II 3mL/200mg	
5 x HR-XC, HR-XA, HR-XCW, HR-XAW	<b>730723</b> (45 μm)	<b>730726</b> (85 µm)	30

Further sizes and adsorbent weights for all HR-Xpert phases on request!



All our CHROMABOND® SPE-phases:

www.mn-net.com

Please contact us ...

Telephone: +49 (0) 2421 969 175 e-mail: tech-chroma@mn-net.com

www.mn-net.com \_\_\_\_\_\_15\_\_\_\_\_



#### More literature needed? Please contact us for:



SPE Info brochure (english)

Catalog (english)



ΜN Chromatography Catalog (german/english/ french)

NUCLEODUR® HPLC Catalog (english)

HPLC Application Guide (english)

Application Guide (english)

... or please visit our online application-database at www.mn-net.com/apps

















#### local distributor

www.mn-net.com

### MACHEREY-NAGEL



and international: Tel.: +49 24 21 969-0 Fax: +49 24 21 969-199 E-mail: info@mn-net.com

Switzerland: MACHEREY-NAGEL AG

Tel.: +41 62 388 55 00 Fax: +41 62 388 55 05 

France: MACHEREY-NAGEL EURL +33 388 68 22 68 +33 388 51 76 88 Tel.: Fax:

USA:

MACHEREY-NAGEL Inc.

Tel.: +1 484 821 0984 Fax: +1 484 821 1272 E-mail: sales-us@mn-net.com

