


NeurosTM Syringes

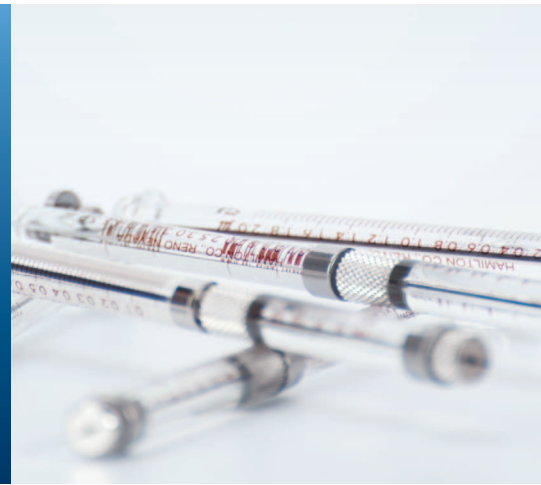

Maximum Control, Minimum Damage


Hamilton NeurosTM syringe technology provides unprecedented functionality for controlled animal injections. An industry first, the Neuros accurately dispenses 50 nL to 100 μ L of liquid through an ultrafine needle protected by a custom-designed sleeve. Developed specifically for the neurosciences, the Neuros minimizes injection site damage by increasing needle rigidity and delivering a microvolume of liquid to an exact location.

Flexible Options for a Variety of Uses


Neuros technology is offered on several models from the 700 MicroliterTM, 1700 Gastight[®] and 7000 Modified Microliter series. These syringes are crafted to have very little to no dead volume which preserves the maximum amount of buffer and solution. Neuros syringes feature a blunt point (point style 3) so the sample is delivered directly to the site; they also feature two different needle sleeves. The sleeve with a blind stop is perfect for cannulated applications and ensures targeted administration with an adjustable penetration depth. The version without a blind stop works best with stereotaxic holders. Both types provide an adjustable needle exposure of 0 to 20 mm.

Benefits of the Neuros Design:

- ▶ Needle rigidity improves insertion path accuracy
- ▶ Minimal tissue damage reduces injection variability
- ▶ Reduced sample loss saves money and materials
- ▶ Fine gauge needle creates smaller injection sites
- ▶ Compatibility with most infusion pumps and stereotaxic holders means an easy integration into existing processes

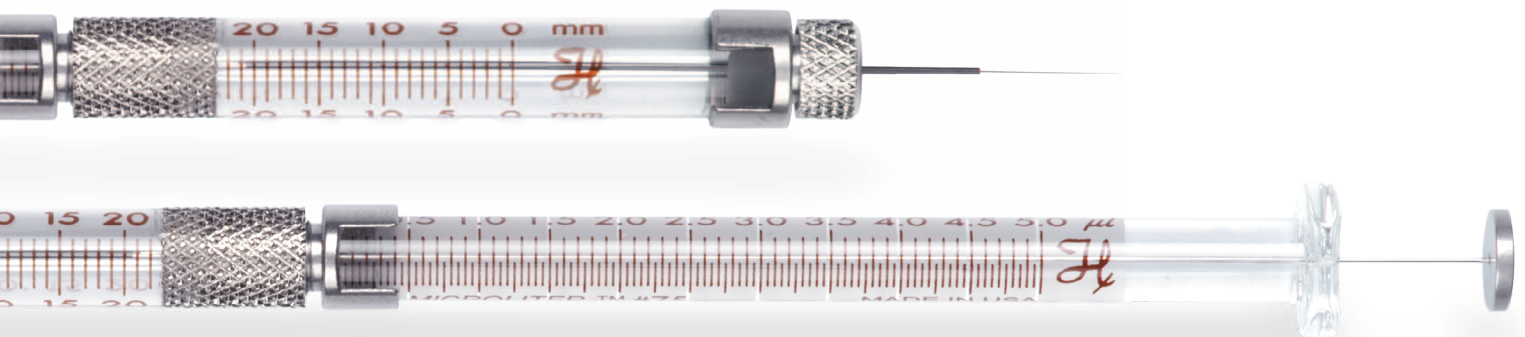

Neuros™ Syringes

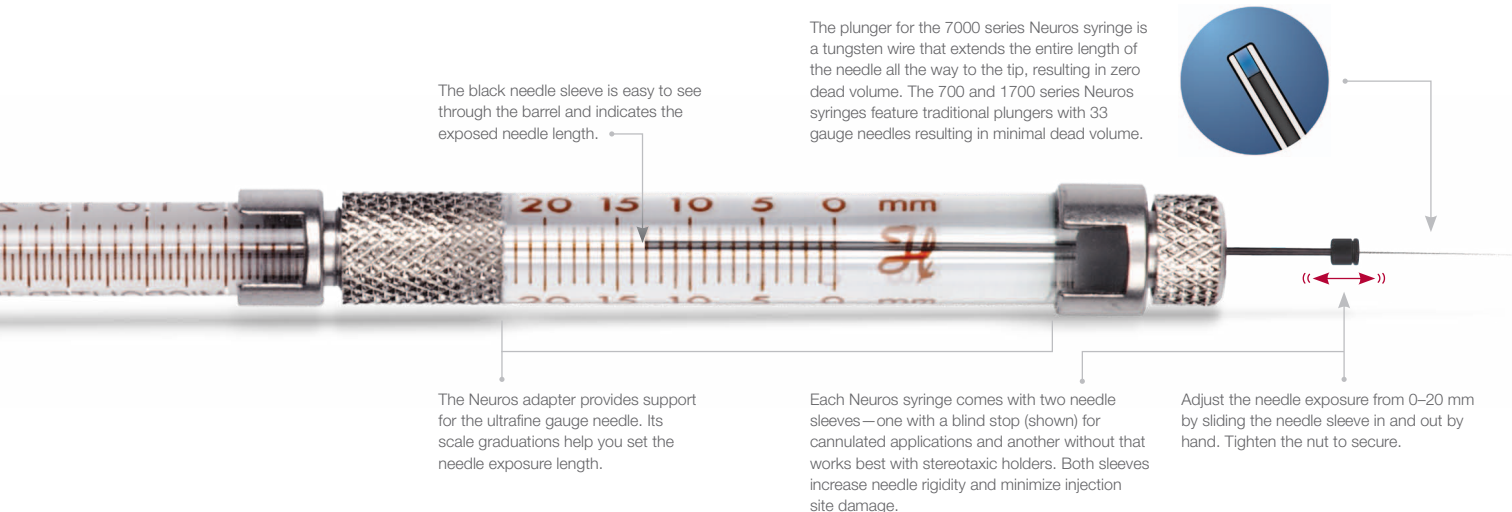

Ordering Information:

Hamilton Part No.	Volume	Description	Gauge	Syringe Type
65457-01	0.5 µL	Model 7000.5 KH Neuros Syringe	32	Modified Microliter
65458-01	1.0 µL	Model 7001 KH Neuros Syringe	32	Modified Microliter
65459-01	2.0 µL	Model 7002 KH Neuros Syringe	30	Modified Microliter
65460-02	5.0 µL	Model 75 RN Neuros Syringe	33	Microliter
65460-05	10 µL	Model 1701 RN Neuros Syringe	33	Gastight
65460-10	25 µL	Model 1702 RN Neuros Syringe	33	Gastight
65460-15	50 µL	Model 1705 RN Neuros Syringe	33	Gastight
65460-20	100 µL	Model 1710 RN Neuros Syringe	33	Gastight

For more detailed information on Neuros technology or to order, visit www.hamiltoncompany.com/neuros.

© 2013 Hamilton Company. All rights reserved.
All trademarks are owned and/or registered by Hamilton Company in the U.S. and/or other countries.
Lit. No. L20076 Rev E

HAMILTON®

Web: www.hamiltoncompany.com

USA: 800-648-5950

Europe: +41-81-660-60-60

Hamilton Americas & Pacific Rim

Hamilton Company Inc.
4970 Energy Way
Reno, Nevada 89502 USA
Tel: +1-775-858-3000
Fax: +1-775-856-7259
sales@hamiltoncompany.com

Hamilton Europe, Asia, & Africa

Hamilton Bonaduz A.G.
Via Crusch 8
CH-7402 Bonaduz, Switzerland
Tel: +41-81-660-60-60
Fax: +41-81-660-60-70
contact@hamilton.ch

To find a representative in your area, please visit hamiltoncompany.com/contacts.